

# incubators

## Percival model I-36LL

### applications

- Designed for applications such as drosophila, culturing, rearing and holding of test organisms as well as bioassays, seed germination and product stability testing
- Many other applications exist for this product  
*Please compare your own requirements to the specifications listed below.*

### percival's intellus ultra controller

- Controls temperature, lighting, humidity (optional) and CO<sub>2</sub> (optional)
- Single-board electronic solid-state design includes 10 key membrane keypad with LED indicators and vacuum fluorescent display
- Programs can be configured to run in real time or elapsed time
- Ramping and non-ramping program methods available for each programming mode
- Multiple programs can be linked creating complex environmental profiles
- Optional Intellus Web Server allows monitoring and controlling of chamber via web browser (requires Internet Explorer 6.0+) (this option allows for remote monitoring and programming of chamber including alerts and current condition updates for up to five e-mail addresses)

*Please refer to [www.percival-scientific.com](http://www.percival-scientific.com) for additional information regarding the control system.*

### lighting system

- Two 20W fluorescent lamps horizontally mounted in pairs above each shelf
- Intensity programmable up to 65  $\mu\text{moles/m}^2/\text{s}$  of light irradiance measured @ 6" from lamps on 2 on/off light events
- Two levels of programming fluorescent lighting
- Programming and control of the lighting is done via Intellus real time controller



### cabinet construction

- Interior constructed of 22-gauge electro-zinc plated steel
- Exterior constructed of 18-gauge exterior electro-zinc plated steel
- Welded seams and joints on outer and inner shells
- Inner shell supported by non-compressing/non-thermal conducting material locking inner liner in place without a metal-to-metal bond to outer case
- Chamber is completely self-contained
- Overall wall thickness is 2" (5.1 cm)
- One 1 1/4" diameter access port on R.H. wall
- Chamber floor equipped with floor drain and hose assembly
- Contains caster assembly and adjustable leveling legs to compensate for floor unevenness in the lab

### insulation

- Woodless construction using CFC free insulation (overall wall thickness is 2" [5.1 cm], ample insulation for maintenance of stated temperature range)

## I-36LL specifications (subject to change without notice)

Temp Range with all lights on	Interior Space volume		Total Shelving Floor Area		Maximum Growing Height		Exterior Dimensions						Light Intensity 6" from lamps unless otherwise noted	Tiers
	°C	ft <sup>3</sup>	m <sup>3</sup>	ft <sup>2</sup>	m <sup>2</sup>	in	cm	width		depth		height		
4-44±0.5	29.7	0.8	21.6	2	10	25.4	33.5	85.1	33.6	85.4	77.2	196.1	65	4

# incubators Percival model I-36LL

## door

- One door opening 29.2" x 57.5" (74 cm x 146 cm) provides full access to the chamber interior (magnetic gasket provides a tight seal to door frame)

## interior space

- 29.7 ft<sup>3</sup> (0.8 m<sup>3</sup>) with work area of 21.6 ft<sup>2</sup> (2.0 m<sup>2</sup>) provided on four tiers

## shelving

- Four tiers of white epoxy coated steel wire shelving (each shelf is 27"D x 28.8"W [68.6 cm x 73 cm])
- Shelves are supported by shelf clips allowing ½" vertical adjustments
- Maximum clearance between shelves is 10" (25.4 cm) per tier with all four shelves installed

## finish

- Interior and exterior painted with highly reflective, environmentally friendly, high temperature baked white powder coating

## refrigeration

- ¼ h.p. self-contained air-cooled condensing unit with hot gas bypass system for continuous compressor operation, extended life and close temperature control (this continuous running condensing unit ensures precise temperature control by alternately cycling refrigerant and hot gas to coil; this also prolongs life of compressor, and eliminates risk of ice build up in coil)
- Top mounted refrigeration system allows easy access for maintenance (e.g. cleaning)
- As heat is rejected, it rises and is dissipated into room without having any effect on inside temperature of cabinet
- Solenoid valves have extended stem for quiet and long life operation
- Ceiling mounted evaporator coil incorporates twin air circulation fans in aluminum housing (heat rejection to ambient [standard chamber] = 3800 BTU/hr.)

## temperature range

- 4°-44°C (±0.5°C) lights on and 2°-44°C (±0.5°C) lights off

## temperature safety limit controls

- (Experiment Protection) Adjustable high and low temperature controls, audible alarms, and visual indicators provided
- Controls shut down all power to the chamber, activating alarms (when the temperature returns to the normal range the system will automatically reset)

## humidity control (optional)

This section outlines the H1 option

- Additive humidity control of higher than ambient to 80% (±10%) lights on for set temperatures between 15° to 30°C

## humidity control (optional, continued)

- Humidity control of higher than ambient to 90% (±10%) lights off for temperatures between 15° to 30°C
- Extended humidity ranges available  
See other specification sheets or consult factory for additional information. If humidity system is selected as an option, a de-mineralized water supply is required which terminates to a ½" MPT connector.

## options (most popular)

- Phenolic Coated Coils (required for drosophila research) (Q9)
- Glass door (Q12)
- Intellus Ultra Web Server (C9)
- Communications Software (C9+)
- Intellus Ultra with Touchscreen and Internet capabilities (C10)
- Pan-type humidifier with Electronic RH Sensor (H1)
- Pan-type humidifier and dehumidifier with Electronic RH sensor (H3)
- Ultrasonic Humidifier with advanced RH Sensor (H11)
- Dehumidification via independent dehumidifying coil with reheat heaters and Ultrasonic Humidifier (H12)
- Ultrasonic Humidifier with Electronic RH sensor (H14)
- CO<sub>2</sub> enrichment package
- Door with observation window and cover (Q2)
- Door with fresh air ports (Q1)
- Self-contained water-cooled condensing unit
- Dry alarm contacts
- Dimmable lighting (closed loop with PAR light sensor) (Q22)
- Dimmable lighting (open loop control) (Q23)
- Extended temperature ranges available  
See other catalog sheets or consult factory for additional accessories.

## convenience receptacles

- Two 115/1/60 convenience receptacles provided inside chamber

## electrical service requirements

- 115/1/60 - 12 amps (total) for standard chamber (power cord and grounded plug provided)



Donauwörther Str. 9  
86637 Wertingen  
Tel.: +49 (0) 8272 6430 60  
Fax: +49 (0) 8272 6430 61

info@plantclimatics.de  
www.plantclimatics.de